
Bedienungsanleitung • Mode d'emploi • Instruction

# OptiMIC

MIKROSKOPE  
MICROSCOPES

**MA 113 / 116 / 118 / 119 / 121 / 130**

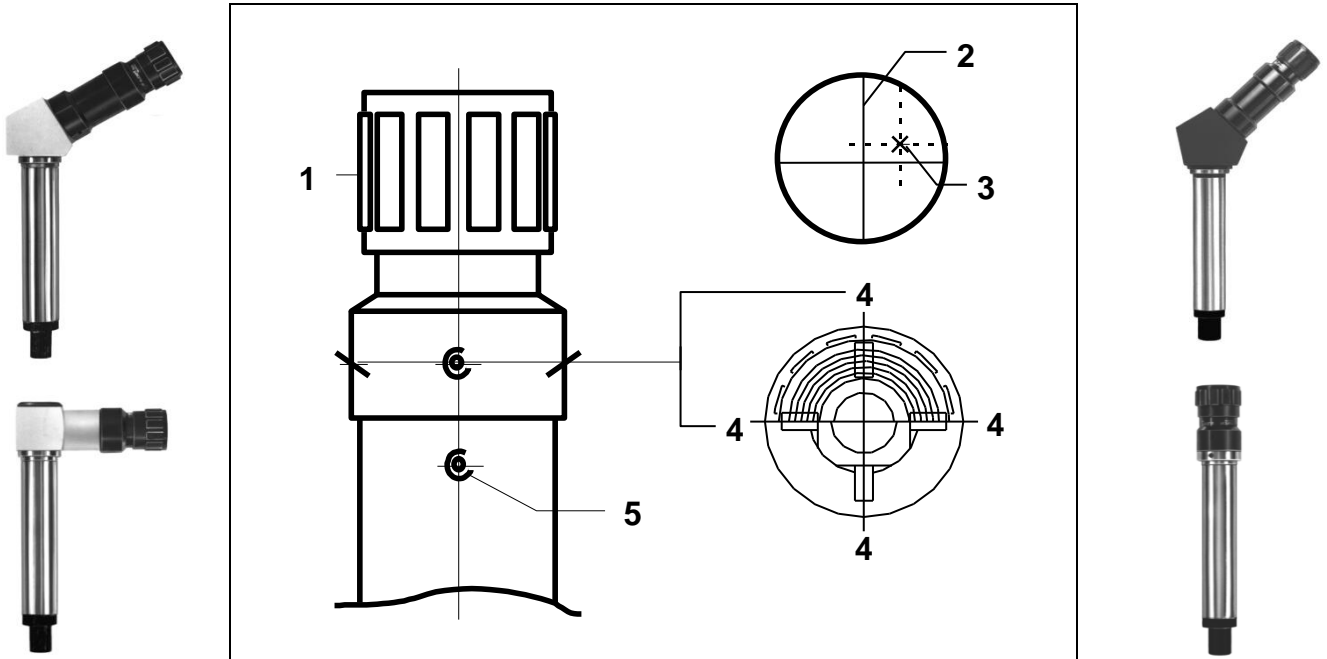## MIKROSKOPE MA 113 / 116 / 118 / 119 / 121 / 130

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---

## *The microscope is a precision instrument and must be treated very carefully.*

- 1) Unpack carefully microscope and accessories. If necessary wipe optic with a lint free clean cloth.
- 2) **Mounting :**
  - Put the microscope in its stand.
  - Turn the diopter ring **(1)**, till the graticule is net for your eye.
  - Adjust the sharpness of the image by a reference piece or a centering point (move the microscope, the reference piece, or the centering point up or down).
- 3) **Eventual centering of the microscope:**

The microscope has been carefully centered and checked before dispatch. It is recommended however to re-check its centering before use as vibration during transport may have caused a misalignment of optical elements.  
If a recentering is necessary, see point 4 „centering of the graticule“.
- 4) **Centering of the graticule:**

The eventual centering fault, i. e. distance between the crosswire **(2)** and the controlpiece **(3)** (reference piece or centering point), can be corrected with the crosswire **(2)** and in moving the 2 screws **(4)** horizontal or vertical.  
**(Attention: The 2 screws (4), horizontal or vertical, has to be moved together, i. e., loose one of the screws and on the same time tight the screw opposite).**

### **Orientation of the graticule:**

Loose the screw **(5)**. Turn the ocular-head **(1)** and after the adjustment tight the screw **(5)**. Re-check the centering of the graticule and if necessary, repeat point 4 „centering of the graticule“.

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